Instructions to collect hair samples

Hair samples can be taken by anyone who has received adequate prior training, and has the materials required.

Materials required:

✓ Card stock with 1 cm on grid

✓ Painters tape, ¼" wide

✓ Ziploc bag / envelope

✓ Surgical clamp (mosquito)

- ✓ Hair grips/clip
- ✓ Comb


✓ Scissors

✓ Gloves (optional, but recommended)



Step 1

• Make a horizontal parting in the hair using the clips, at the back of the head (posterior vertex).

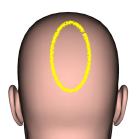



Step 2

 Separate a section of hair (thickness of a pencil, 0.5cm) about 4 cm above the skull base.



Step 3

- Clamp a lock of hair 2cm from the root, if the length permits.
- If the hair is too short or not very thick, you can make a zigzag movement through the hair before tightening the hair with the clamp, to avoid leaving a visible hole on the scalp.
- You can also take hair from several locations in the same region in order to obtain the same amount of hair.

Step 4

- In the clamp, the width of hair strand should be 1cm wide. Measure with the help of the grid on the card.
- Using the scissors, cut the hair as close to the scalp as possible.






Step 5

- While the hair is still in the clamp, place the section of hair on the card with the root upwards.
- Using painter's tape, secure the lock of hair on the card.
- Repeat as needed if the hair is collected in several places.
- It is important that the root of the hair is placed according to the arrows on the card.

Step 6

- Identify the sample with a label placed on card.
- Place the card in a Ziploc bag, making dure that the sample is dry.
- Seal the bag and put a label on it.

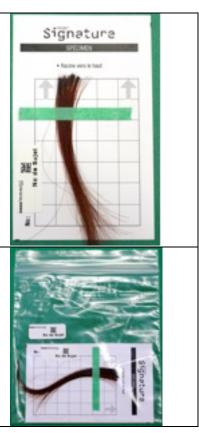

Store samples at room temperature, in a dry and dark place. Sample bags should not be placed tightly together. A card filling cabinet is ideal.

Make note of the following informations for each participant: Hair coloring and/or treatment, frequency of washing, season of collection, topical and oral medication, mental and physical health.